



**USE & CARE GUIDE**

**MED2 Series  
REFRIGERATORS & FREEZERS  
FOR MEDICAL & LABORATORY STORAGE**

If you need service or call with a question, please have this information ready. We advise you to keep this *Use & Care Guide* in your possession.

Model: \_\_\_\_\_ Serial number: \_\_\_\_\_

Options added: \_\_\_\_\_

Date of purchase: \_\_\_\_\_ Dealer: \_\_\_\_\_

NOTE: The alarm/thermometer included in this unit has a NIST calibrated probe that requires re-calibration every two years. The NIST calibration expiration date is included on the device. In order to maintain your NIST certification, you will need to re-calibrate your probe at that time.



# MODELS

## All-Refrigerators

- **FFAR24LMED2**  
*Compact all-refrigerator with automatic defrost operation*
- **FF28LWHMED2**  
*Countertop all-refrigerator with automatic defrost operation*
- **FF511LBIMED2 / FF511LBIMED2ADA**  
*Undercounter all-refrigerator with automatic defrost operation*
- **FF511LBI7MED2 / FF511LBI7MED2ADA**  
*Commercially listed undercounter all-refrigerator with automatic defrost operation*
- **FF7LBIMED2 / FF7LBIMED2ADA**  
*Undercounter all-refrigerator with automatic defrost operation*
- **FFAR10MED2**  
*Mid-sized all-refrigerator with automatic defrost operation*
- **SCR600LBIMED2 / SCR600LBIMED2ADA**  
*Undercounter glass door all-refrigerator with automatic defrost operation*

## All-Freezers:

- **FS24LMED2**  
*Compact all-freezer with manual defrost operation*
- **FS24L7MED2**  
*Commercially listed compact all-freezer with manual defrost operation*
- **FS30LMED2**  
*Countertop all-freezer with manual defrost operation*
- **FS30L7MED2**  
*Commercially listed countertop all-freezer with manual defrost operation*
- **FS407LBIMED2 / FS407LBIMED2ADA**  
*Undercounter all-freezer with manual defrost operation*
- **FS407LBI7MED2 / FS407LBI7MED2ADA**  
*Commercially listed undercounter all-freezer with manual defrost operation*
- **VT65MLBIMED2 / VT65MLBIMED2ADA**  
*Undercounter all-freezer with manual defrost operation*
- **VT65MLBI7MED2 / VT65MLBI7MED2ADA**  
*Commercially listed undercounter all-freezer with manual defrost operation*

# Table of Contents

<b>Model Numbers</b> .....	<b>1</b>
<b>Appliance Safety</b> .....	<b>3</b>
<i>Proper disposal of an old appliance</i> .....	3
<b>Installation Instructions</b> .....	<b>4</b>
<i>Key Eject lock</i> .....	4
<b>Location Requirements</b> .....	<b>5 – 6</b>
<i>Location</i> .....	5
<i>Recommended grounding method</i> .....	5
<i>Leveling your appliance</i> .....	6
<i>When using the appliance</i> .....	6
<i>Energy-saving tips</i> .....	6
<b>Appliance Care</b> .....	<b>7 – 8</b>
<i>Defrosting</i> .....	7
<i>Cleaning your appliance</i> .....	7
<i>Changing the light bulb</i> .....	8
<i>Loss of power</i> .....	8
<i>Moving the appliance</i> .....	8
<b>Temperature Selection and Monitoring</b> .....	<b>9 – 11</b>
<i>Digital thermostat</i> .....	9
<i>Thermometer/alarm</i> .....	10 –12
<b>Keypad Lock</b> .....	<b>13</b>
<b>Operation of Specific Models</b> .....	<b>14</b>
<b>Troubleshooting</b> .....	<b>15</b>
<b>Limited Warranty</b> .....	<b>16</b>

# Appliance Safety

Your safety and the safety of others are very important to us.

We have provided many important safety messages in this manual and on your appliance. Always read and obey all safety messages.

You can be seriously injured or even killed if you don't follow instructions.

Safety messages will tell you what the potential hazard is, how to reduce the chance of injury, and what can happen if the instructions are not followed.



## IMPORTANT SAFETY INSTRUCTIONS

**WARNING: To reduce the risk of fire, electric shock or injury when using your appliance, follow these basic precautions:**

- Plug into a grounded 3-prong outlet.
- Do not remove ground prong, nor use a 3/2 adapter.
- Do not use an extension cord.
- Be sure not to plug into outlets controlled by wall switches.
- Disconnect power before servicing.
- After servicing, replace all parts and panels before operating.
- Use nonflammable cleaners.
- Keep flammable materials and vapors, such as gasoline, away from appliances.
- Use two or more people to move and install appliance.

### **Proper disposal of an old appliance:**

**DANGER! Risk of child entrapment!**

**Old appliances can represent a suffocation hazard to children.**

- Remove doors from your old appliance.
- Leave shelves in place so children cannot easily climb inside.

If an old appliance contains refrigerants, it must be recycled by a licensed service or disposal company.

# Installation Instructions



## **Excessive Weight Hazard !**

**Use two or more people to move and install your appliance.**

**Failure to do so can result in back or other injury.**

**ALLOW APPLIANCE TO SIT UPRIGHT FOR 24 HOURS BEFORE TURNING ON, TO ALLOW THE COMPRESSOR LUBRICANT TO DRAIN.**

1. Remove packaging materials.
2. Remove tape and glue residue from surfaces before turning on the appliance. Rub a small amount of liquid dish soap over the adhesive with your fingers. Rinse with warm water and dry.
3. Do not use sharp instruments, rubbing alcohol, flammable fluids, or abrasive cleaners to remove tape or glue. These products can damage the surface of your appliance.
4. When Moving Your Appliance:  
Your appliance is heavy. When moving the appliance for cleaning or service, be sure to protect the floor. Always pull the appliance straight out when moving it. Do not wiggle or “walk” the appliance when trying to move it, because floor damage can occur.
5. Clean Before Using:  
After you remove all of the packaging materials, clean the inside of your appliance before using it. See cleaning instructions in *Appliance Care*.
6. Some upright models have leveling legs. Level as necessary for best operation.
7. If your appliance has casters, after moving into place, lock the brakes on those casters supplied with brakes. Always unlock before moving the appliance for cleaning.

### **Key Eject lock (some models):**

To lock or unlock the appliance, you must hold the key in the lock while turning. The lock is spring loaded, so the key cannot be left in the lock.

If the key is lost, contact your dealer or SUMMIT APPLIANCE for a replacement. Always give the model and serial number of your appliance when requesting a replacement key. Keys for most models can be ordered online at [www.summitappliance.com/support](http://www.summitappliance.com/support)

### **A NOTE ABOUT DOOR SWING:**

Due to their wiring and controls, the doors on all MED2 units **cannot be reversed in the field**. When ordering in the future, always specify whether you need a right hand or left hand (option “LHD”) door swing. All units ship with a right hand door swing unless specified differently at the time of order.

## Location Requirements



**Keep flammable materials and vapors, such as gasoline, away from appliances. Failure to do so can result in fire, explosion or death.**

### Location:

**IMPORTANT:** Do not install the appliance near an oven, radiator, or other heat source, nor in a location where the temperature will fall below 40°F (5°C). Do not place in direct sunlight.

Check the strength of the floor before installing your appliance. It should hold the weight of a fully loaded appliance. Allow at least 4" (10 cm) of airspace at the top and on both sides of the appliance and also 4" between the back of the appliance and the wall. If the appliance is going to be placed next to a wall, leave extra space on the hinge side so the door can open wider than its width.

NOTE: Allow 4" (10 cm) clearance at the top and sides and a ¼" (0.64 cm) clearance between appliances. **IF YOUR MODEL HAS THE LETTERS "BI" IN THE SUFFIX, IT IS DESIGNED FOR BUILT-IN USE AND REQUIRES ONLY MINIMAL CLEARANCES EXCEPT AT THE BACK WHERE AT LEAST 2" OF CLEARANCE SHOULD BE ALLOWED.**



### Electrical Shock Hazard

- **Plug into a grounded 3-prong outlet.**
- **Do not remove ground prong.**
- **Do not use an adapter.**
- **Do not use an extension cord.**

**Failure to follow these instructions can result in fire, electrical shock or death.**

### Recommended grounding method:

A 115 Volt (AC only) 60 Hz grounded electrical supply with a 15- or 20- amp fuse or circuit breaker is required. It is recommended that a separate circuit serving only your appliance be provided. Use outlets that cannot be turned off by a switch or pull chain. Do not use an extension cord.

## **Leveling your appliance (some models):**

Locate the two front leveling legs that may be taped among the packing materials. (In many models these legs are already mounted into the underside of the appliance.)

1. Screw these legs into the front holes on the bottom of the appliance. The back of the appliance rests on two fixed supports. (Built-in models also have rear level legs that can provide more precise leveling).
2. Adjust the front legs just enough to lift the cabinet up off the floor.  
**Note:** If your model is equipped with front rollers, the front legs should be lowered enough to lift the rollers off the floor to ensure the appliance does not roll forward when the door is opened.
3. Place a leveling tool on top of the appliance, first side to side, then front to back. Turn leg counter-clockwise to raise a corner. Turn leg clockwise to lower a corner.
4. Repeat Step 3 until appliance is level.
5. After appliance is fully loaded, recheck with level as in Step 3.
6. If your appliance is equipped with casters, the two front casters have brakes to prevent rolling. Keep brakes locked, and unlock for moving only.

## **When using the appliance:**

In order to ensure proper temperature, you need to permit proper airflow in the appliance. Do not block off any air vents. If the vents are blocked, airflow will be hindered, and temperature and moisture problems may occur. Always position contents to allow some airflow between shelves or baskets.

Allow the appliance to cool at least 6 to 8 hours before adding contents. Allow 24 hours for temperature to stabilize.

NOTE: When the appliance is first started, the compressor will run constantly until the appliance is cooled. It may take 6 to 8 hours (or longer), depending on room temperature and the number of times the appliance door is opened. After the appliance has cooled, the compressor will turn on and off as needed to keep the appliance at the proper temperature.

**For optimum temperature performance, it is recommended that this unit be at least 30% full during use. If you are only using this unit to store small loads, we recommend adding some water bottles or freezer packs to ensure temperature stability.**

The outside of your appliance may feel warm. This is normal. The appliance's design and main function is to remove heat from packages and air space inside the appliance. The heat is transferred to the room air, making the outside of the appliance feel warm. *(In some models the condenser that gets warm is just inside the side walls, making these walls warm to the touch.)*

## **ENERGY-SAVING TIPS**

- The more airflow provided around an appliance, the less energy it will use.
- If your appliance requires manual defrosting, keep ice build-up to less than 1/2".
- Keep doors open a minimum of time, and shut securely.

# Appliance Care

## Defrosting:

Refrigerator models listed in this manual defrost automatically. However, freezer models require manual defrosting. Before defrosting the unit, remove the contents of the freezer, then set the thermostat to OFF (the compressor will stop working). Leave the door open until ice and frost are completely melted. To speed up defrosting, you can place a container of warm water (about 125°F) into the cabinet. Soak up the melt water with clean toweling or a sponge and be sure the interior is dry before turning the thermostat back to its normal setting.

- NOTE: It is not advisable to heat the interior of the freezer directly with hot water or a hair dryer while defrosting since this can deform the cabinet interior.

**NEVER USE A METAL OBJECT TO REMOVE ICE FROM THE REAR WALL. IF NEEDED, USE A PLASTIC ICE-SCRAPER.**

## Cleaning the appliance:

**NOTE:** To prevent the buildup of odors, clean the appliance about once a month. Follow the procedure below for cleaning:

1. Disconnect from power.
2. Remove any parts that can be taken out of the cabinet, such as drawers.
3. Using a clean sponge or soft cloth, hand wash, rinse and dry removable parts and interior surfaces thoroughly. Use a mild detergent in warm water. DO NOT use abrasive or harsh cleansers such as window sprays, scouring cleansers, flammable fluids, cleaning waxes, concentrated detergents, bleaches or cleansers containing petroleum products. DO NOT use harsh cleaning tools such as paper towels or scouring pads, as these can scratch or damage the surfaces of the refrigerator.
4. For stainless steel and painted metal exterior surfaces, use a clean sponge or soft cloth and a mild detergent in warm water. DO NOT use abrasive or harsh cleaners, or cleaners containing chlorine, as these can scratch or damage the material. Using a soft cloth, dry thoroughly.
5. To help with removing odors, you can use a mixture of warm water and baking soda (2 tbs to 1 qt of water) to wash the interior walls.
6. After cleaning, replace all of the parts that were removed from the appliance.
7. The condenser coils, located on the back or behind the base grille, should also be cleaned regularly. This may be as often as every other month. Cleaning the condenser coils can help reduce how much energy the appliance uses. When the condenser coils are dusty or dirty, use a vacuum cleaner with an extension attachment.
8. After cleaning the appliance, plug in or reconnect power.

***NOTE: For CSS Models (completely stainless steel), only use cleaners especially made for stainless steel.***

### **Changing the light bulb (some models only):**

NOTE: Not all appliance bulbs will fit your appliance. Be sure to replace the bulb with one of the same size, shape, base size and wattage. You can purchase bulbs from [www.summitappliance.com/support](http://www.summitappliance.com/support)

1. Unplug appliance or disconnect power.
2. Remove the light shield. Most models do not require tools.
3. Replace the light bulb with an appliance bulb of the same wattage. Replacement LED bulbs are available from our parts department.
4. Replace the light shield.
5. Plug in appliance or reconnect power.

### **Loss of power:**

In the event of power loss, unplug the appliance. Once power is restored, allow 30 minutes for the refrigeration system to equalize. During this time, keep the door closed as much as possible.

**Remember:** A full appliance stays cold longer than a partially filled one.

### **Moving the appliance:**

1. Remove stored items and tape shelves, drawers and trays in place.
2. Unplug the appliance.
3. Clean it thoroughly.
4. Tape the appliance shut and tape the electrical cord to the cabinet.

When you get the appliance to its new location, refer to the *Installation Instructions* and *Temperature Selection* sections for information on installation and setting the controls.

## Temperature Selection & Monitoring



*Note: details may vary from image*

### Digital Thermostat

#### **Temperature Parameters:**

**MED2 all-refrigerators** come with digital thermostats that are preset at 4°C, with a range of 2° to 8°C.

**VT65MLBIMED2 and VT65MLBIMED2ADA** all-freezers are preset to -20°, with a range of -25° to -10°C.

All other **MED2 all-freezers** are preset to -18°C, with a range of -20° to -10°C.

If you find it necessary to change the temperature setting for your digital thermostat, please refer to the following instructions.

#### **How to Adjust the Temperature:**

##### **To see the SET POINT:**


The default display for your digital thermostat is the internal cabinet temperature. To view the unit's set point, push and immediately release the *SET* key. The display will flash, showing the *set point* value. If no other actions are taken in 6 seconds, the display will revert to the internal air temperature.

##### **To change the SET POINT:**

While the *set point* value is blinking, press the UP or DOWN arrow to change the set point to your desired temperature. Press the *SET* button again to exit the adjustment. The display will revert to the internal air temperature. **PLEASE NOTE:** The temperature displayed on the digital thermostat is the internal air temperature. Your alarm/thermometer (installed on the left in the control panel) provides a temperature readout based on the buffered probe in order to better simulate the temperature of what is actually stored inside the unit (as opposed to the air temperature itself). As a result, these readings may not always display the same temperature reading.

**NOTE: The minimum and maximum range of your thermostat is preset and locked by us, so it cannot easily be accidentally moved outside of the recommended temperature ranges.**

## Thermometer/Alarm

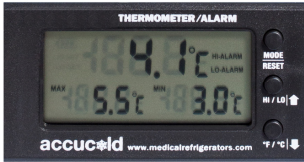

*Note: Details may vary from image*

Every MED2 unit is equipped with a NIST calibrated Thermometer/Alarm that simultaneously displays minimum, maximum, and current interior temperatures and provides a visual and audible signal when the temperature inside the unit falls outside the set temperature range.

**NOTE ABOUT CALIBRATION:** All MED2 units ship with a two-year **certificate of calibration** in the alarm/thermometer box that comes inside your unit. Keep this certificate stored safely for your records. Accucold offers re-calibration services through their ISO 17025:2005 certified calibration laboratory. It is recommended to have your probe re-calibrated every two years (or one year, depending on your state's health department guidelines). For more information on how to have your device re-calibrated, contact us at [calibration@summitappliance.com](mailto:calibration@summitappliance.com).

The temperature sensor, or probe, is enclosed in a glycol-filled bottle that is stored inside your unit in a special shelving compartment.

In some units, the center shelf is fixed in place and cannot be removed (see image). This is because the glycol-filled bottle that contains the temperature sensors for your alarm and thermometer is best stored in this location. The center of your refrigerator or freezer will provide the most accurate temperature readings. For proper temperature monitoring, keep the bottles in place and do not remove this shelf.



## **Thermometer/alarm instructions (cont'd):**

### **Function of Controls:**

**MODE/RESET Key:** Press ONCE to reset Max/Min temperature record. When the unit is sounding an alarm, press ONCE to turn off the beeping sound.

**HI/LO Key:** Press ONCE to display high alarm setting. Press TWICE to display low alarm setting. Press THREE TIMES to display current temperature reading.

**°F/°C Key:** Press this button to change the temperature display scale from Fahrenheit (°F) to Celsius (°C).

**Battery Installation:** Remove battery compartment cover located on the rear of the unit and remove insulation strip from battery compartment. Following the polarity (+/-) diagram in the battery compartment, install the battery. Replace the battery cover. The unit will beep and all segments of the LCD will be activated.

### **Programming the Alarms:**

The thermometer displays the internal temperature of the unit being monitored, as well as the maximum and minimum temperatures reached. The maximum and minimum readings reflect the highest and lowest temperatures since the battery was inserted or since the MODE/RESET key was last pressed. If the temperature rises above or falls below the set temperature range, the alarm will sound. By default, the high alarm temperature setting is 8°C (46.4°F) and the low alarm temperature setting is 2°C (35.6°F). To reset these limits, follow these instructions:

**High Alarm Temperature Setting:** Press the HI/LO key ONCE, then press MODE/RESET to enter the **High Alarm** temperature setting mode. Use the Up and Down arrows to adjust the temperature accordingly. Press MODE/RESET once again to confirm the setting.

**Low Alarm Temperature Setting:** Press the HI/LO key TWICE, then press MODE/RESET to enter the **Low Alarm** temperature setting mode. Use the Up and Down arrows to adjust the temperature accordingly. Press MODE/RESET ONCE to confirm the setting.

By default, the low and high alarms should be audible. To turn alarms on/off:

**Audible High Alarm Setting:** To make sure the high alarm is audible, press HI/LO key ONCE (to activate HI-SET on the screen). Then press the ALERT ON/OFF key ONCE to activate the HIGH-ALARM beeping sound. Press again to deactivate.

**Audible Low Alarm Setting:** To make sure that the low temperature alarm is audible, press the HI/LO key TWICE to activate LOW-SET on the screen. Then press the ALERT ON/OFF key ONCE to activate the LOW-ALARM beeping sound. Press again to deactivate.

NOTE: Set high and low alarm temperature limits above and below the refrigerator/freezer operating temperature to prevent false alarms. For operating temperature ranges, see *Temperature Parameters* on page 9.


NOTE: For an activated high alarm and low alarm buzzer, the display should indicate HI-ALARM and LO-ALARM (see illustration on previous page).


**More details about the features and usage of the Thermometer/Alarm can be found in the User Manual accompanying the unit.**

## Keypad Lock (optional on most models)


**Factory-set default password:** The factory-set default password is 0123. If the default



password is lost, press the  key (located inside the battery box) for 30 seconds, then wait for the warning sound which signifies the factory-set default password has been restored.

**To set the password:** While door is open, press the  key (located inside the battery box) until it jingles, enter the 0123 password, then # within 10 seconds to confirm. Enter your password (1-16 digits), then press # again to confirm. A total of 10 passwords can be set.


NOTE: Each password starts with a number from 0 to 9 in order. The first password starting with 0 is the default password. For example, use 1xxx, then 2xxx etc. for 10 users.

**Changing password:** While the door is open, press the  key (located inside the battery box) until it jingles, enter the old password and press # within 10 seconds to confirm. Enter the new password, then press # to confirm.

NOTE: The first digit cannot be changed.

**To open the door:** While the door is closed, press the START/END key, enter the password, then press START/END again to confirm.

**To lock door manually:** Press #.

**Function setting of automatic lock:** Setting of the auto-lock time: Press the  key (inside the battery box), enter a number (any number between 10-99), then press # to confirm. The auto-lock time can be set between 10-99 seconds after opening the door. If the time is set less than 10 seconds, the door will not lock automatically, only manually.

**Deleting the door-opening passwords:** Press the START/END key, enter default password followed by 0000, then press # and all the door-opening passwords will be deleted except for the default password.

NOTE: To access the battery box you need to remove only the two small screws on the top and bottom of that cover. Do not remove the 4 screws holding the keypad and cover to the appliance door.

## Operation of Specific Models

**All-Refrigerators: FF24LMED2, FF28LWHMED2, FF511LBIMED2, FF511LBI7MED2, FF511LBIMED2ADA, FF511LBI7MED2ADA, FF7LBIMED2, FF7LBIMED2ADA, FFAR10MED2, SCR600LBIMED2, SCR600LBIMED2ADA**

These models are all-refrigerators (no freezer compartment) with automatic defrost. All have cold plates on the rear wall, and the coldest spots in the refrigerator will be in the rear and toward the bottom. All have interior fans to promote even cooling. In the rear bottom above the step, there is a drain line to allow defrost water to exit the refrigerator. Keep this drain line clear, or icing inside the refrigerator will occur. FF7, FF511L, FFAR10, and SCR series models have lights that are not door-activated, and are controlled either by a rocker switch on the top right or a light switch inside. The light inside these units will stay on until switched off manually. While LED lighting will not generate heat, it is not recommended to leave the light on for long periods of time.

***NOTE:** Models with BI in the model number are designed for built-in operation and need very little clearance on the sides and above the units. All other models require at least 4 inches of airflow on the sides and above the units. All models, including built-in units, operate best when 2 inches is provided between the rear wall and the back of the unit.*

**All-Freezers: FS24LMED2, FS24L7MED2, FS30LMED2, FS30L7MED2, FS407LBIMED2, FS407LBI7MED2, FS407LBIMED2ADA, FS407LBI7MED2ADA, VT65MLBIMED2, VT65MLBIMED2ADA, VT65MLBI7MED2, VT65MLBI7MED2ADA**

These are manual defrost all-freezers. No internal fans are present, and ice will build up along the cooling coils. Periodically defrost the ice (recommended when the ice level reaches a thickness of 1/4 to 1/2 inch). **NEVER USE A METAL INSTRUMENT** as you can damage the cooling coils, which will void your warranty and often cannot be repaired. Before defrosting the unit, remove the contents of the freezer, then set the thermostat to OFF (the compressor will stop working) and disconnect the power. Leave the door open until ice and frost are completely melted. To speed up defrosting, you can place a container of warm water (about 125°F) into the cabinet. Soak up the melt water with clean toweling or a sponge and be sure the interior is dry before turning the thermostat back to its normal setting.

*For built-in (BI) models and clearance information, see NOTE above.*

# Troubleshooting

Try these solutions first to avoid the cost of an unnecessary service call.

***PROBLEM:*** *The appliance is not operating.*

- Is the power supply cord unplugged? Plug into a grounded 3-prong outlet.
- Has a fuse blown or a circuit breaker been tripped? Replace the fuse or reset the circuit.
- Is the thermostat incorrectly set? See *How to Adjust the Temperature* on page 9.

***PROBLEM:*** *The motor is constantly running, or running too much*

- Is it hotter than usual where the appliance is located? At normal temperatures, the motor will typically run between 40% and 80% of the time. The motor will run even longer when the unit is in an unusually warm environment.
- Have a large number of items been added to the appliance? The more that is added, the longer it will take for the unit to cool down. During this process, the motor will run longer.
- Is the door frequently opened? Warm air entering the appliance causes the motor to run longer. Keep the door closed as much as possible.
- Is the thermostat set correctly? See *How to Adjust the Temperature* on page 9.
- Is the door completely closed? Push the door firmly shut.
- Are the condenser coils dirty? This prevents proper air transfer, and causes the motor to work harder. Clean the condenser coils. See *Cleaning the Appliance* on page 7, and refer to #7.
- Does the unit have proper clearance? Models with "BI" in their name require 2" of rear clearance and minimal clearance on the sides and top. All other units require at least 4" of clearance on all sides for proper operation.

***PROBLEM:*** *The temperature in the cabinet is too warm*

- Have you just added a large number of items to the appliance? This will warm the interior of the unit until it is able to cool down to the correct temperature.
- Is the thermostat set correctly? See *How to Adjust the Temperature* on page 9.

***PROBLEM:*** *There is excessive moisture inside the appliance*

- Is the door opened frequently? If so, humidity from the room will enter the appliance and cause moisture to build up. This will happen even more if the air in the room is very humid.

***PROBLEM:*** *The door is hard to open*

- Is the door gasket dirty or sticky? Clean the gasket and the surfaces that it touches. Rub a thin coat of paraffin wax on the gasket after cleaning.

# Limited Warranty

## **2-YEAR LIMITED WARRANTY FOR APPLIANCE**

Within the 48 contiguous United States, for two years from the date of purchase, when this appliance is operated and maintained according to instructions attached to or furnished with the product, warrantor will pay for factory-specified parts and repair labor to correct defects in materials or workmanship. Service must be provided by a designated service company. Outside the 48 states, all parts are warranted for two years from manufacturing defects. Plastic parts, shelves and cabinets are warranted to be manufactured to commercially acceptable standards, and are not covered from damage during handling or breakage.

## **5-YEAR COMPRESSOR WARRANTY**

1. The compressor is covered for 5 years.
2. Replacement does not include labor.

### ***ITEMS WARRANTOR WILL NOT PAY FOR:***

1. Service calls to correct the installation of your appliance, to instruct you how to use your appliance, to replace or repair fuses or to correct wiring or plumbing.
2. Service calls to repair or replace appliance light bulbs or broken glass shelves. Consumable parts (such as filters) are excluded from warranty coverage.
3. Damage resulting from accident, alteration, misuse, abuse, fire, flood, acts of God, improper installation, installation not in accordance with electrical or plumbing codes, or use of products not approved by warrantor.
4. Replacement parts or repair labor costs for units operated outside the United States.
5. Repairs to parts or systems resulting from unauthorized modifications made to the appliance.
6. The removal and reinstallation of your appliance if it is installed in an inaccessible location or is not installed in accordance with published installation instructions.

## **DISCLAIMER OF IMPLIED WARRANTIES – LIMITATION OF REMEDIES**

CUSTOMER'S SOLE AND EXCLUSIVE REMEDY UNDER THIS LIMITED WARRANTY SHALL BE PRODUCT REPAIR AS PROVIDED HEREIN. IMPLIED WARRANTIES, INCLUDING WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE, ARE LIMITED TO ONE YEAR. WARRANTOR SHALL NOT BE LIABLE FOR INCIDENTAL OR CONSEQUENTIAL DAMAGES. SOME STATES DO NOT ALLOW THE EXCLUSION OR LIMITATION OF INCIDENTAL OR CONSEQUENTIAL DAMAGES, OR LIMITATIONS ON THE DURATION OF IMPLIED WARRANTIES OF MERCHANTABILITY OR FITNESS, SO THESE EXCLUSIONS OR LIMITATIONS MAY NOT APPLY TO YOU. THIS WARRANTY GIVES YOU SPECIFIC LEGAL RIGHTS AND YOU MAY ALSO HAVE OTHER RIGHTS, WHICH VARY FROM STATE TO STATE.

The following warning is required for goods sold in the state of California:



**WARNING: This product can expose you to chemicals including Nickel (Metallic) which is known to the State of California to cause cancer.**

For more information go to [www.P65Warnings.ca.gov](http://www.P65Warnings.ca.gov)

Note: Nickel is a component in all stainless steel and some other metal components.

**accucold**