

# GTR-AS / Adjustable Speed Tube Rocker Quick Start Guide

## What's in the box:

- Quick Start Guide
- Tube Rocker
- Power Cables (1 each: US plug, UK plug, EU plug)
- 10 Position Silicone Tray Insert



**IMPORTANT:** The tube rocker must not be overloaded or damage may occur. Please do not exceed 1kg (2.2 lbs).

## INSTALLATION & OPERATION

### Operating the Unit

- Carefully unpack and place the unit on a stable non-slip surface.
- Allow approximately 4 inches surrounding the unit for safe operation.
- Connect the power adaptor using the correct plug for your power supply.
- Turn power switch on.
- Rotate the speed control knob to the desired speed setting.