

# LAB REFRIGERATORS & FREEZERS

## GENERAL USE / QUESTIONS

---

***Q: What is the weight limit of the shelves?***

A: It is not recommended to exceed 75 Lbs. per shelf. Due to shelf weight distribution, it is not recommended to have more than 8 shelves per door.

***Q: Should I leave my equipment operating when I am not using it?***

A: This unit is designed to operate continuously. Leaving the refrigeration system in operation, (even if not in use) may extend the life of the equipment, and reduce the chance of refrigeration issues that may occur if the unit is not in operation for long periods of time.

***Q: What ambient conditions is this unit designed for?***

A: This unit is designed for:

- Indoor use
- Altitude up to 2000m.
- Ambient Temperatures 15°C to 32°C ( 60°F TO 85°F )
- Recommended humidity range of 30% to 90%.

***Q: What electrical conditions is this unit designed for?***

A: Individual unit specifications may vary. See the documentation on the back of the unit for clarification on electrical requirements.

***Q: Does my unit require a dedicated electrical line?***

A: This unit does require a dedicated electrical line. See the documentation on the back of the unit for clarification on electrical requirements.

***Q: Does the battery power my unit when power is lost?***

A: No, if a battery was included with your unit, it only powers the alarm system, and will not keep the unit cold during a power outage.

***Q: How often does the battery need to be replaced?***

A: If a battery was included with your unit, it is recommended to replace the battery at least once every 3 years.