


Pipette Shaker

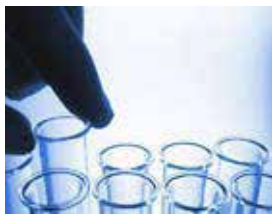

Unpacking and Setup

Carefully unpack the pipette shaker, saving packing materials in the unlikely event you need to return the rocker. Immediately call LW Scientific if you detect any damage. If you need to return the item, please use the original packaging.

Setup

Place Pipette Shaker on a flat, stable surface. Ensure the controls are set in the "OFF" position. Plug into the appropriate electrical outlet.

Controls

- Three position Power Switch: Used to turn the Pipette Shaker on and off. Positions are ON, OFF, and TIMER. ON setting runs continuously.
- Red Power light: Signals the unit is on.

Operation

1. Place samples on the top of the holder
2. Turn on the Power switch to ON or TIMER position

Specifications

Speed:	2500 RPM
Power:	110V / 60Hz (220V available)
Height:	158.75mm (6.25 in)
Width:	133.35mm (5.25 in)
Weight:	1.9Kg (4.2 lbs.)

Notes for Safe Use:

- Do not operate unit in or near water or other liquids.
- Unplug unit before cleaning or repairing.
- Do not use if electrical cord is damaged.
- Replace cord if damaged.